



**WATERJACKET CO2
INCUBATORS**

WITH MICROPROCESSOR CONTROL

MODELS: 3507, 3517, 3514, 3524

INSTALLATION AND OPERATION MANUAL

Sheldon Manufacturing Inc. P.O. Box 627 Cornelius, Oregon 97113
EMAIL: tech@Shellab.com **INTERNET:** <http://www.Shellab.com/~Shellab>
1-800-322-4897 (503) 640-3000 FAX (503) 640-1366

TABLE OF CONTENTS

SECTION 1.0	RECEIVING AND INSPECTION
SECTION 2.0	GRAPHIC SYMBOLS
SECTION 3.0	INSTALLATION
SECTION 4.0	CONTROL PANEL OVERVIEW
SECTION 5.0	OPERATION
SECTION 6.0	DATA LOGGER FUNCTION
SECTION 7.0	FYRITE CO2 CHECKING
SECTION 8.0	HEPA CO2 FILTER
SECTION 9.0	MAINTENANCE
SECTION 10.0	TROUBLESHOOTING
SECTION 11.0	PARTS LIST
	UNIT SPECIFICATIONS
	SCHEMATICS

REV 03/06
4861607

These units are CO2 Water Jacketed incubators for professional, industrial or educational use where the preparation or testing of materials is done at approximately atmospheric pressure and no flammable, volatile or combustible materials are being heated. These units are not intended for hazardous or household locations or use.

RECEIVING AND INSPECTION




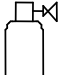







IMPORTANT: READ THIS INSTRUCTION MANUAL IMMEDIATELY!

Your satisfaction and safety require a complete understanding of this unit, including its proper function and operational characteristics. Be sure operators are given adequate training before attempting to put the unit in service. NOTE: This equipment must be used only for its intended application; any alterations or modifications will void your warranty.

- 1.1 Inspection:** The carrier, when accepting shipment, also accepts responsibility for safe delivery and is liable for loss or damage claims. On delivery, inspect for visible exterior damage, note and describe on the freight bill any damage found and enter your claim on the form supplied by the carrier.
- 1.2** Inspect for concealed loss or damage on the unit itself, both interior and exterior. If any, the carrier will arrange for official inspection to substantiate your claim. Save the shipping crate until you are sure the unit has been delivered in good condition.
- 1.3 Return Shipment:** If for any reason you must return the unit, contact your customer service representative for authorization and supply nameplate data. For information on where to contact Customer Service please see the manual cover.
- 1.4 Accessories:** Make sure that all of the equipment indicated on the packing slip is included with the unit. Carefully check all packaging before discarding. Each chamber comes equipped with 3 shelves, 4 shelf standards, 1 humidity pan and a power cord. All units are shipped with 4 leveling feet and supply hose kit.

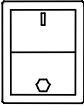GRAPHIC SYMBOLS

Your incubator has been provided with a display of graphic symbols which is designed to help in identifying the use and function of the available user adjustable components.

- | | | |
|------|---|--|
| 2.1 |  | Indicates that you should consult your manual for further description or discussion of a control or user item. |
| 2.2 |  | Indicates " Temperature ". |
| 2.3 |  | Indicates " Over-temperature safety ". |
| 2.4 | °C | Indicates " Degrees Centigrade ". |
| 2.5 | CO ₂ | Indicates " Carbon Dioxide ". |
| 2.6 |  | Indicates " Gas " (CO ₂ for this unit). |
| 2.7 |  | Indicates " AC Power ". |
| 2.8 |  | Indicates " Water Jacket Low ". |
| 2.9 |  | Indicates " ON/I " and " OFF/O ". |
| 2.10 |  | Indicates " Relative Humidity ". |
| 2.11 | %RH | Indicates " Percent Relative Humidity ". |
| 2.12 |  | Indicates " Potential Shock Hazard " behind this protective partition. |
| 2.13 |  | Indicates " Protective Earth Ground ". |
| 2.14 |  | Indicates " Unit should be recycled " (Not disposed of in land-fill) |

INSTALLATION

Local city, county, or other ordinances may govern the use of this equipment. If you have any questions about local requirements, please contact the appropriate local agency. Installation may be performed by the end user.

Under normal circumstances this unit is intended for use indoors, at room temperatures between 5° and 40°C, at no greater than 80% Relative Humidity (at 25°) and with a supply voltage that does not vary by more than 10%. Customer service should be contacted for operating conditions outside of these limits.

CAUTION: Make sure that the incubator is located in its intended position and level before filling the water jacket. Directions for filling the jacket are in Section 5.2.

- 3.1 Power Source:** The power supply must be properly grounded (earthed) and correctly sized to match the unit nameplate rating. **The supply voltage must match the nameplate voltage within +/- 10%.** These units are intended for 50/60 Hz application. If supplied with a detachable cord set, plug the female end into the inlet on the unit and the male plug into the supply. Assure that units requiring a fuse have a fuse installed. This fuse may be at the inlet or a part of the cord set male plug.
- 3.2 Location:** When selecting a site for the unit, consider conditions which may affect performance, such as heat from steam radiators, ovens, autoclaves, etc. Avoid direct sun, fast-moving air currents, heating/cooling ducts, and high-traffic areas. To ensure air circulation around the unit, allow a minimum of 10cm of clearance between the incubator and surrounding walls, or obstructions to free air flow.
- 3.3 Lifting / Handling:** These units are heavy and care should be taken to use appropriate lifting devices that are sufficiently rated for these loads. Units should only be lifted from their bottom surfaces. Doors, handles and knobs are not adequate for lifting or stabilization. The unit should be completely restrained from tipping during lifting or transport. All moving parts, such as shelves and trays should be removed and doors need to be positively locked in the closed position during transfer to prevent shifting and damage.
- 3.4 Leveling:** The unit must sit level and solidly. Leveling feet are supplied and must be installed in the four holes in the bottom corners of the unit. With the feet installed and the unit standing upright, each foot can be raised by turning it in a counterclockwise direction. Adjust the foot at each corner until the unit stands level and solid without rocking. If the unit must be moved, drain all of the water from the unit and turn the leveling feet in (clockwise) all the way to prevent damage while moving.
- 3.5 Cleaning:** The unit chamber should be cleaned and disinfected prior to use.

Your operating conditions and appropriate protocol will determine the correct procedure for decontamination. A typical decontamination procedure that is adequate for many situations has been described below. As well, certain steps are listed that will help reduce the likelihood of contamination and the necessity of decontamination. Whatever process is appropriate, it needs to be done on a regularly scheduled basis. Depending on usage and protocol, this may be monthly, quarterly or otherwise. Regardless of the decontamination procedure used, certain precautions will need to be taken:

- A. Always disconnect the unit from the electrical service when cleaning. Assure all volatile or flammable cleaners are evaporated and dry before reconnecting the unit to the power supply.
- B. Special care should be taken when cleaning around sensing heads to prevent damage.
- C. If you use chlorine-based bleaches or abrasive cleaners this will modify the stainless steel interior finish. DO NOT USE hard tools such as metal wire brushes or steel wool. Use non-abrasive cleaners and soft tools such as plastic brushes.

TYPICAL DECONTAMINATION PROCEDURE

1. Remove the humidity pan every week and autoclave, or wash with soap and water then disinfect with 70% alcohol solution. Replace in the incubator with fresh, DISTILLED water.
2. Flush the sample port tubing with 70% alcohol solution. Replace any lines that have a green color.
3. Remove the door gaskets, clean and disinfect. Clean and disinfect all mounting grooves for the door gaskets.
4. Remove all shelves, shelf supports, shelf standards and shields. Autoclave, or wash and disinfect as described in item 1.
5. Wash and disinfect all interior surfaces.
6. Give special attention to cleaning and disinfecting all access ports, air bleeds, shaft holes, electrical pass-through and any other passages into the chamber.
7. Replace all air and CO₂ filters every six months or when noticeably dirty on the upstream side. CO₂ filters are located in the shadow box just behind the GAS IN fitting and in line with the CO₂ tubing kit. Units with flow meters require operator- supplied air pumps. These pumps should be checked for proper air filtering.


OPERATION FOR MINIMIZING CONTAMINATION

1. Keep the outside of the incubator, including the air in the laboratory, as clean as possible. This is particularly important for units placed directly on the floor. Do not place incubators near doors, air vents or other areas of high air movement or traffic.

2. The floor around the unit needs to be clean. Units that are placed on the floor should be mounted higher – typically on a caster platform – for ease of moving the unit during cleaning and access to the back of the unit.
3. Minimize the number of times access is made to the chamber during normal operation.
4. Do not depend on the use of antibiotics to maintain uncontaminated conditions, as this is an inadequate technique for sterilization. Preferably use aseptic techniques as described above for maintaining sterile conditions in the incubator.

3.6 Shelves and Interior Parts: Shelving, shelf supports and a humidity pan are supplied with each unit. **See Figure 1**, for proper placement of these accessories.

Figure 1

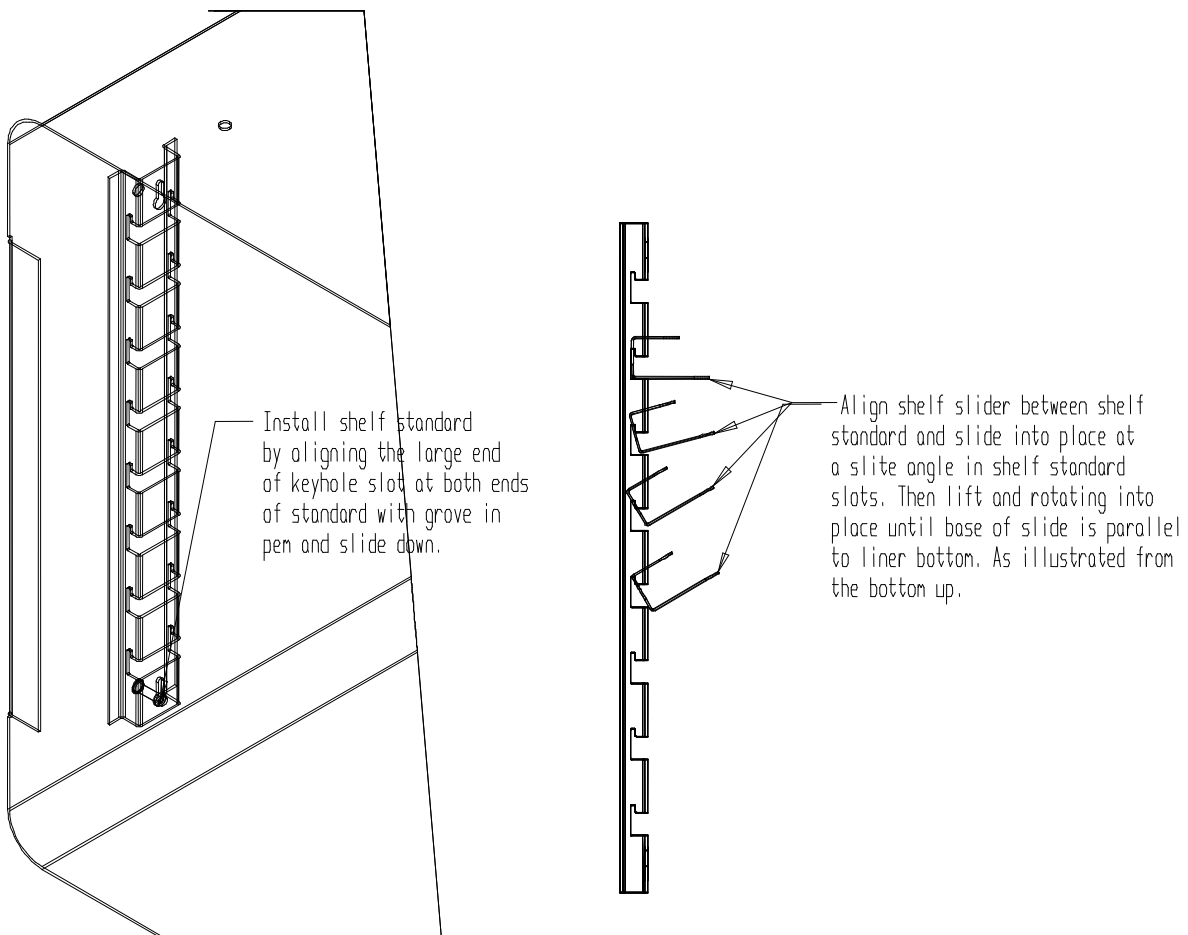

CONTROL PANEL OVERVIEW (See Figure 2)

All controls are located on the front panel. Units with a detachable cord have a fused inlet located at the top rear of the control panel. This inlet has a recessed male plug, fuse and an EMI filtering system designed to filter out electrical interference. This inlet also prevents any internally generated interference from feeding out to the power grid.

- 4.1 Power Switch:** The I/O (ON/OFF) switch controls all of the power for the incubator and must be in the I/ON position before any systems are operational. Both temperature and CO₂ displays will illuminate when the power switch is in the ON position.
- 4.2 Water Low Light:** This pilot lamp is on whenever the internal water jacket float switch has been tripped to the closed position. When the water drops low enough the float switch closes the circuit turning on the indicator light.
- 4.3 Over Temperature Light:** This pilot lamp is on whenever the Over Temperature Safety thermostat has been activated and taken control of the element. During normal operating conditions this indicator should never be on.
- 4.4 Main Temperature Control:** This controller is marked in C° and indicates the actual temperature within the chamber to .1°C. The UP/DOWN buttons are used for inputting the set point, calibrating the display, and muting or unmuting the audible alarm. The HIGH and LOW alarm indicators will light whenever there is an alarm condition associated with the temperature within the chamber. The MUTE indicator will light whenever the audible alarm has been deactivated. Some alarm conditions have a delay programmed in.
- 4.5 Heating Light:** This pilot lamp is on whenever the Temperature Controller is activating the heating element.
- 4.6 CO₂ Control On Water Jacket Incubators:** The control sets the rate and the limit of injection of CO₂ into the incubator based on input value and some variables. The input value is set by the operator. This is set at 5% CO₂ in production. The variables for units with TC sensors are the amount of humidity in the chamber, and the amount of time the control has been trying to inject CO₂.


At start-up, the incubator must be warmed up to operating temperature. This is set in production at 37°C. During warm up, the water tray for humidity should be filled and placed inside. This allows the humidity to reach 95% before CO₂ is injected. The TC type CO₂ sensor requires 95% humidity to correctly measure

the CO2 levels.

The CO2 should be turned OFF during warm up. Failure to do so will cause the control overshoot the set value input by the operator, the injection value can be reset by setting the CO2 levels down until CL shows on the display. Then resetting the CO2 control to the desired setting resets the injection rate and limit values to that which was input by the operator.

- 4.7 **Injection Light:** This pilot lamp is on whenever the CO₂ controller is injecting CO₂ into the chamber.
- 4.8 **Humidity Display:** This digital display is marked %RH and indicates the relative humidity in the chamber.
- 4.9 **Datalog Communications Interface:** This controller provides six (6) parameters to externally log data of all the front panel displays. The MODE button is used to select the parameter, and the UP/DOWN buttons are used to change the parameter's value.
- 4.10 **Over Temperature Safety Control (OTP):** Located front left of the control panel, this is a hydraulic thermostat that is wired between the Main Temperature Controller and the heating element and functions as an override control. If at any time the Main Temperature Control fails in the ON position and the temperature in the chamber rises above its set point, the OTP is activated and maintains temperature at its own set point. Note that the HEATING indicator will continue to function under the control of the OTP.
- 4.11 **Interface Ports:** Located on the front left of the control panel opposite the Over Temperature Safety, these three (3) ports are where the logged data can be externally accessed through serial and parallel functions.
- 4.12 **CO₂ Sample Port:** This is located on the front left side of control panel. A sample can be drawn to measure the CO₂ content in the chamber at this port.

Figure 2

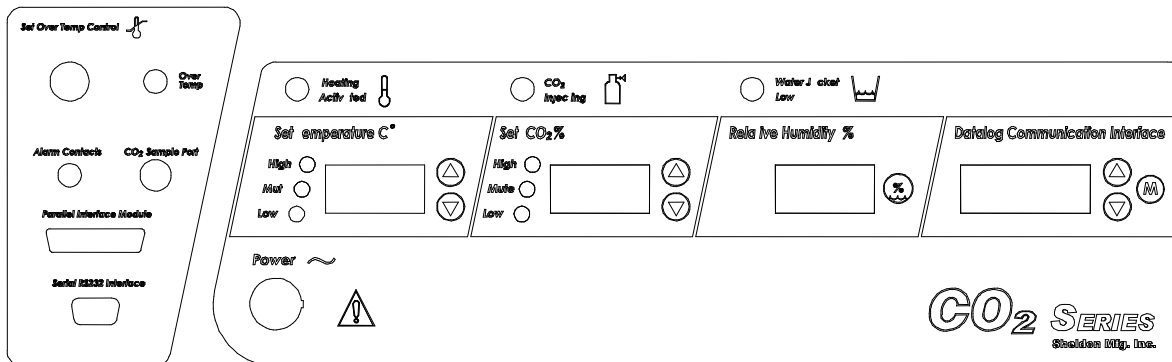

OPERATION

READ THIS SECTION IN ITS ENTIRETY BEFORE ATTEMPTING OPERATION!

IMPORTANT: *In order to ensure the proper operation of this incubator, it is essential that you regularly verify the accuracy of the temperature and CO₂ controls. Temperature can be confirmed by use of a thermometer inside of the chamber. CO₂ can be confirmed with the use of a Fyrite Gas Analyzer or other equipment such as a gas chromatograph. These must be checked at least on a weekly basis. Any questions should be directed to a technical service representative.*

- 5.1 Power up the unit by attaching the cordset (supplied with the accessories) to the unit inlet and the building power supply. Attach the female end of the cordset into the fused and recessed male plug on the rear of the unit and the male plug end of the cordset into the building supply receptacle. Turn the power switch to the I/ON position.
- 5.2 **Filling and Draining:** The front panel of the incubator has a WATER LOW warning light. Fill the water jacket with water by means of the fill port located at the top right rear of the unit marked FILL WJ HERE. Stop filling when the WATER LOW warning light goes off. When full, the unit holds approximately 19 U.S. gallons (72 liters). To drain the unit, hook a suitable hose to the siphon tube marked SIPHON HERE located to the left of the fill port. Start a siphon and allow unit to empty into a suitable service drain or catchment area.

WATER RECOMMENDATIONS FOR USE IN WATER JACKET INCUBATORS

The addition of a dis-similar metal anode in Water Jacket incubators will allow the use of good-quality tap water to fill the Jacket. The anode is located in a threaded port near the fill port on the rear of the unit. The anode will dissolve as it is in contact with mineral salts or dissolved gasses in the tap water. The anode should be removed and checked for erosion every year. When the anode appears to be over 50% eroded, a new one should be purchased from Sheldon Mfg., Inc. (Replacement Part #0260500)

Tap water with a hardness of more than 30 parts per million (PPM) or 1.5 grains per gallon (GPG) will require assessment of the anode on a yearly basis. Dissolved gasses should be no more than 120 PPM or 7 GPG. The Ph of the water should test to between 6.0 and 8.5.

When in doubt, it is advisable to use DISTILLED WATER ONLY.

5.3 Setting the Controls: Refer to the Control Panel Overview Section 4.6 of this manual.

5.4 Calibrating the Controls:

- A. Push and hold both the UP and DOWN buttons simultaneously until the control displays CP (for temperature) or CO (for CO₂) after the display will start to blink.
- B. While the display is blinking the UP and DOWN buttons can be used to adjust the display to match the actual condition in the incubator chamber.
- C. If no buttons are pressed within five (5) seconds the blinking will stop and the display will revert to showing the process or actual parameter within the incubator chamber.

5.5 Humidification: Humidification of the unit is achieved by evaporation of water from the chamber humidity pan placed in the bottom of the incubator. By filling this stainless steel reservoir pan with **DISTILLED WATER ONLY** and allowing this water supply to heat and evaporate, near saturation humidity is obtained. Do not use plastic, glass or other metals. Only 300 series stainless metals are acceptable for this reservoir pan. Do not use corrosive chemicals including copper sulphate or chlorine in the pan or chamber as damage may occur. **Use Distilled Water Only. DO NOT USE DEIONIZED WATER!**

Use of disinfecting chemicals in the pan can change the surface tension of the reservoir water thus preventing evaporation and proper humidification of the chamber. Water in the pan should be changed and the pan cleaned at least once a week to help control contamination and to maintain proper surface tension.

5.6 Temperature Monitoring: To insure that the incubator is operating at the desired temperature, an accurate temperature indicator such as a certified reference thermometer should be placed in the incubator chamber. Try to place the thermometer in the center of the chamber and spaced off the shelf (taping the thermometer down to a petri dish cover is a method that raises it off the shelf and prevents it from rolling and hiding the scale from view). For ease of reading the thermometer would be better if they were electronic with a remote display so they could be read without opening the doors.

5.7 Initial Control Settings: Set controls to the set point you want your incubator to operate at.

- A. Set the OTP to its maximum position clockwise. After making initial control settings allow the incubator to warm up and stabilize for 24 hours.


- B. All units are shipped set for 37°C and CO₂ OFF. If other conditions are desired follow the instructions in section 5.3 for resetting the controls.

5.8 Temperature Calibration: It is important that the calibration be performed in the following order: Main temperature, and then the Over Temperature Safety.

- A. During main temperature calibration it is important that the inner glass door not be opened for any reason, and that the outer heated door be opened as little as possible.
- B. After the incubator has been running for 24 hours and the temperature display is stable at the set point, read the temperature in the chamber from the reference thermometer, and if necessary, recalibrate the display using the instructions in section 5.4.
- C. After the main temperature display has restabilized and maintained set point for several hours, check the actual temperature again. If it does not match the display repeat the calibration.
- D. After the Main Temperatures are set and adjusted, the Overtemperature Safety Control needs to be set. Do this by turning the OTP Control Knob counterclockwise until the Overtemperature Light comes on (this will blink in sequence with Heating Activated Light). Next, slowly turn the knob clockwise until the light goes off, then turn the knob another 8 degrees approximately. This should set the Overtemperature 1 degree above the Main Temperature set point.

5.9 CO₂ Supply System and Control System: The CO₂ system is rated for pressures between 10 to 20 PSI which should never be exceeded at any time. Prior to any attempt to hook up or run the CO₂ control system, the water jacket must be filled and the temperatures adjusted as noted in section 5.7.

Figure 3

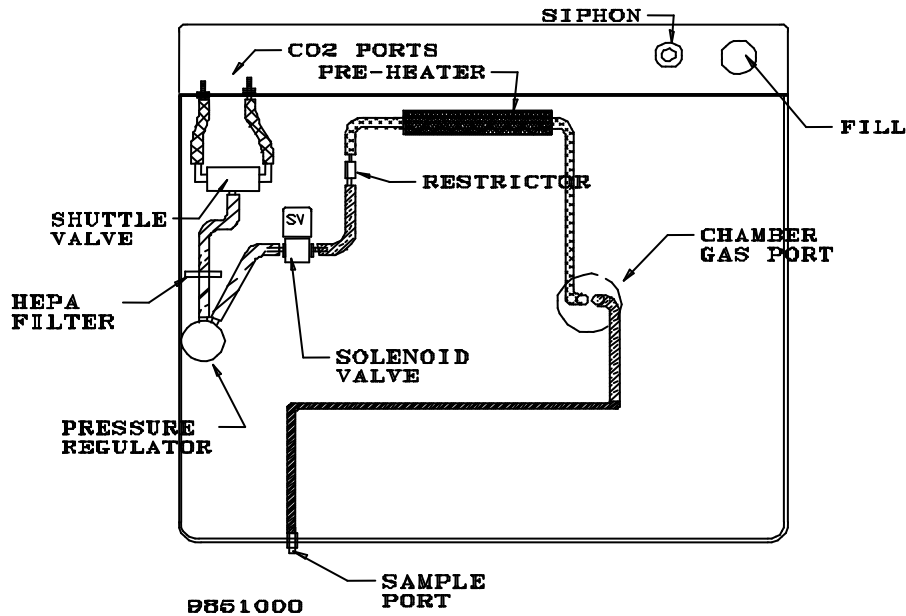

- A. The CO₂ inlet fittings are located at the back of the unit near the top (see **figure three**). They are marked CO₂ TO CHAMBER. You have the option of using one or two tanks. A supply hose with in-line CO₂ filter (supplied with your accessories) connects from each fitting to a CO₂ tank and regulator. If you plan to use only one tank, it does not matter which fitting you connect the supply hose to, and the remaining fitting is left open. If you plan to use two tanks, set the output pressure of one tank to 15 PSI and the other tank to 20 PSI. The internal tank switch will use up the 20 PSI first and then switch over to the 15 PSI tank. It does not matter which tank is attached to which fitting. When the 20 PSI tank is empty its output pressure gauge will indicate 15 PSI. At this time the empty tank can be replaced. Note that there are no alarms to indicate when the tanks are empty, so operators should monitor gas consumption.
- B. It is highly recommended that a good quality 0-60 PSI output range **DUAL STAGE** pressure regulator be used on the CO₂ tank. The dual stage regulator will have two pressure gauges. The high pressure gauge (0-4000 PSI) will indicate the pressure within the tank. The low pressure gauge (0-60 PSI) will indicate the output pressure on the supply hose to the incubator. Do not use single stage regulators. They will not supply a stable pressure. Just because a regulator has two gauges does not mean that it is a dual stage regulator. **Insure that you are using a dual stage regulator.**

- C.** It is normal for the high pressure gauge on your regulator to start out reading 800 to 1000 PSI with a full tank. The reading will drop to 500 to 800 PSI quickly and will stay there for most of the duration of the tank. At the end of use the pressure will drop quickly to zero to indicate that the tank is completely empty. Pure CO₂ is in a liquid state in the tank, and a constant vapor pressure is generated in the tank above the liquid level. The CO₂ is drawn off of the top as a gas. The same vapor pressure is maintained as long as any liquid is left in the tank. When the last of the liquid has evaporated into gas then the pressure will drop rapidly as the gas is drawn off.
- D.** Only medical grade CO₂ should be used in your incubator. The use of any lesser grade may damage your incubator and void your warranty.
- E.** The micro processor CO₂ control system interprets the information from the CO₂ sensor, displays the CO₂ concentration directly on the digital display, reads the CO₂ set point and controls the percentage of CO₂ in the incubator chamber.
- F.** There are two types of CO₂ sensors available, thermal conductivity (TC) and infrared (IR). The 3507 series operates with TC sensors, and the 3517 series operates with IR sensors.
- G.** The TC controlled unit is referred to as an indirect sensor and operates under the principle that CO₂ is 1.5 times denser than air. If you measure the changes in density they can be interpreted as changes in CO₂ content. CO₂ Gas, therefore, is not directly measured, hence indirect. This method of measurement is accurate, reliable and stable as long as all factors that influence air density are maintained unchanging. Factors that effect density are humidity and temperature. When the door of the incubator is opened, the humidity is reduced, the temperature is lowered and the TC control system loses stability until the conditions in the incubator chamber have recovered to original conditions.
- H.** The IR controlled unit is known as a direct sensor and operates under the principal that a certain frequency of infrared light is absorbed by CO₂. The more CO₂ is present the more light is absorbed. Its accuracy is consistent no matter what the conditions are in the incubator. The IR controller utilizes an infared spectrophotometer to directly read the flow of CO₂ gas into the chamber.
- I.** It is recommended that the accuracy of your CO₂ control system be monitored by measuring the actual CO₂ concentration on a weekly basis with a Fyrite or other measuring device. This should be done when the chamber has not been disturbed for an extended period of time, i.e. after the weekend, or first thing on Monday morning.

5.10 Setting the CO₂ Control: Attach the supply hose to the incubator and turn on the CO₂ supply. Set the CO₂ control to the desired set point using the procedure described in section 5.3.

5.11 Adjusting CO₂ Display: After the incubator has had several hours to stabilize at CO₂ set point, measure the actual CO₂% with a Fyrite. If there is any difference between the Fyrite and the display, use the procedure described in section 5.4 for making the display match the Fyrite. See Section 7.0 for Fyrite use.

NOTE: When using the Fyrite, insure that gas is not being injected while the reading is being taken. Changing the CO₂ set point to 0.0 prior to taking the sample and change the set point back to the desired value after the use of the Fyrite is finished.

5.12 Audible Alarms: The temperature control and the CO₂ control are both equipped with visual and audible alarms.

A. Temperature Alarms: The temperature control has alarm indicators for high and low conditions that are activated whenever the actual temperature is 1° above or below the set point.

B. CO₂ Alarms: The CO₂ controller has alarm indicators for high and low alarm conditions that are activated whenever the actual CO₂% is 1% above or below the set point.

C. Audible Alarms: Both controls are equipped with audible alarms that are activated when either of the HIGH or LOW indicators are activated. The audible alarms can be muted for a single alarm occurrence by pressing and holding down either the UP or DOWN button for several seconds until the alarm mutes. There is a built in delay of approximately 15 minutes on the occurrence of a LOW alarm. This time delay prevents the audible alarm from activating every time the door is opened and the temperature and CO₂ drops.

DATA LOGGER FUNCTION

The data logger provides ways to externally indicate or log all the incubator front panel displays at intervals ranging from 1 to 99 minutes. The logger contains a real time clock/calendar to date and time stamp each logged reading.

The logger indicates and logs all front panel display readings in single lines of text. This data is logged when the present logging interval has expired or one of the monitored displays has changed its alarm state. Logged data has a time and date field that corresponds with the loggers real time clock.

This logged data can be externally accessed through ports provided on the front of the incubator for serial and parallel functions, as well as a BNC contact closure plug for extended alarm condition(s). The parallel port is compatible with any Epson or IBM style printer. The serial port operates at 9600 baud with 8 data bits, 1 (one) stop bit and no parity and connects with a computer through the (DB-9 DCE) RS232.

- 6.1 Program the Data Logger:** Select the parameter to adjust by pressing and holding the "MODE" button pad until the left two display digits show "LP" indicating the logger has entered the set mode. The parameter that can be changed will blink and its value can be changed by pushing the UP or DOWN arrow pads.
- 6.2** Pressing and holding the MODE pad for approximately one second will cause the logger to step to the next parameter. The table on the following page shows the indications for each parameter. If no buttons are pressed for fifteen (15) seconds, the panel display will revert to the normal operation mode.

**THE CONTROL PANEL DISPLAYS
SIX PARAMETERS, SHOWN BELOW, WHICH CAN BE CHANGED USING THE
CONTROL PANEL BUTTONS**

FRONT PANEL MODE	LEFT TWO DIGITS	RIGHT TWO DIGITS	COLON
NORMAL OPERATION	SHOW HOURS OF CURRENT 24 HR TIME	SHOW MINUTES OF CURRENT TIME	BLINK
SET LOGGING PERIOD	SHOW "LP"	BLINK LOGGING INTERVAL IN MINUTES. CAN BE VERIFIED FROM 0-99 MINUTES .	OFF
SET HOURS	BLINKS HOURS PORTION OF CURRENT TIME. CAN BE VARIED FROM 00 TO 23	ARE BLANK	ON
SET MINUTES	ARE BLANK	BLINK MINUTES PORTION OF CURRENT TIME. CAN BE VARIED FROM 0 TO 59	ON
SET DATE	ARE BLANK	BLINK DAY OF MONTH. CAN BE VARIED FROM 1 TO 31	OFF
SET MONTH	BLINK CURRENT MONTH. CAN BE VARIED FORM 1 TO 12	ARE BLANK	OFF
CURRENT YEAR	"19" OR "20"	BLINK LOWER 2 DIGITS OF YEAR	OFF

FYRITE CO₂ CHECKING


A Bacharach FYRITE CO₂ Gas Analyzer is recommended to measure CO₂ concentrations in the incubator chamber. This test instrument is not supplied with the incubator but is readily available from your dealer. Follow the instructions provided with each Fyrite instrument carefully to insure correct and accurate readings.

- 7.1. Press button on top of Fyrite canister to release CO₂ concentration. Tip canister to the side to ensure all fluid is released from top of canister.
- 7.2. Loosen screw on slide scale and align top of fluid with zero on the scale. Tighten screw.
- 7.3. Connect hose and aspirator bulb to unit being tested. The sample port for connection is located on the control panel.
- 7.4. Place the hose sampling cap directly over the plunger valve on top of canister and depress firmly.
- 7.5. With button depressed, squeeze bulb 27 times. On the last squeeze and with bulb still deflated, release hose from button.
- 7.6. Turn Fyrite canister upside down 3 times, each time allowing all fluid to flow to the opposite end of the canister.
- 7.7. Tip canister to the 45 degree position to ensure all fluid has been released from top of canister.
- 7.8. Read CO₂ concentration in %.

NOTE: Your Fyrite indicator will come with a complete set of detailed instructions which should be followed carefully. The fluid used inside this Fyrite instrument is poisonous and corrosive and must not be taken internally. In the event of a spill or accidental body contact with the Fyrite fluid, follow the instructions on the refill bottle carefully.

HEPA CO2 FILTER

The CO2 Incubator design features a HEPA filtration system with a patented copper housing. (Patent No. 6,333,004) As the chamber air is drawn through the filter system, airborne microbes and isolated particulates are not only removed from the air, but destroyed as well. It is through copper oxidation that the contaminants are actually killed. The HEPA filter is easily replaced without the need of any tools and has an efficiency rating of 99.97% at 0.3 microns.

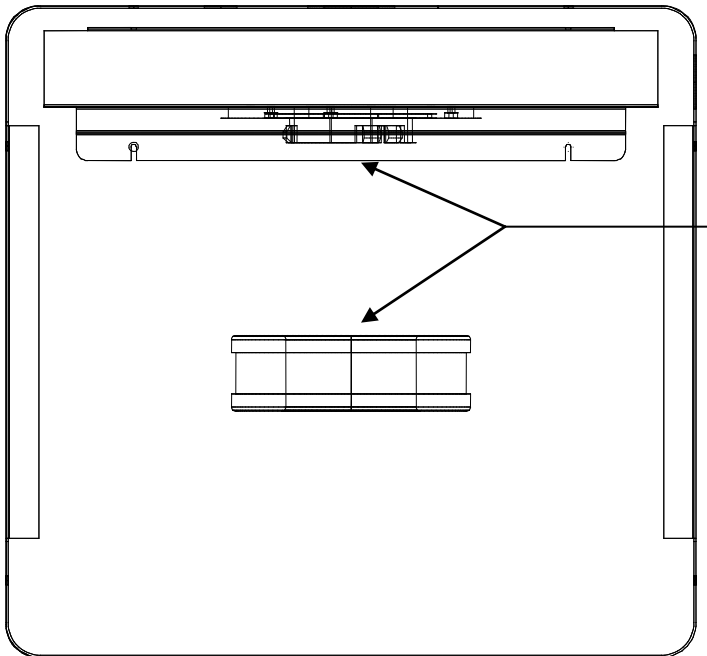

The Hepa filter attaches to a fixed coupling ring with a direct “snap-in” action. Push straight up to the coupling ring to connect. Pull the filter straight down to disengage. Replacement Hepa filters are available from Sheldon Mfg., Inc. Part number 2800517.

MAINTENANCE

NOTE: *Prior to any maintenance or service on this unit, disconnect the service cord from the power supply.*

- 9.1 Cleaning:** Cleaning and decontamination are recommended on a regular basis. To prepare the incubator for cleaning remove all interior parts such as the shelf standards and false top. All stainless steel parts are autoclavable. Please review the Cleaning procedures described in the Installation section for detailed instructions.
- 9.2** Check CO₂ supply periodically; don't let it run out. The internal auto tank switch is not alarmed.
- 9.3** CO₂ supply lines and connections can be checked for leaks with a liquid soap solution. Apply liquid soap solution to all areas of the supply line and observe for bubbles that will indicate a leak.
- 9.4** Keep the CO₂ flow system free of impurities. Erratic CO₂ control is usually traceable to the CO₂ pressure regulator on the tank, impurities in the tank, or impurities in the solenoid valve. Replace the CO₂ in-line filters every six (6) months or when the filters have become noticeably dirty on the upstream side. There is a CO₂ filter attached to the tubing kit, and one inside the shadow box connected to the GAS-IN line.
- 9.5** If the data logger fails to function as specified, check to verify that the port connections are solid and not loose. If the data logger still fails to function properly call Customer Service for assistance.
- 9.6** There is no maintenance required on any controls. If the controls fail to operate as specified, please review Section 9.0 Troubleshooting, prior to contacting customer service.
- 9.7 Storage:** If the incubator is to be turned off for any length of time, dry the chamber and humidity pan thoroughly and leave at room temperature. Failure to do this may cause the interior to become contaminated. No adjustment to controls should be required when restarting the unit. If the unit is to be shut down for transport, drain the water jacket, remove humidity pan and shelves, dry the chamber and disconnect the CO₂ and power supply's. See section 3.3, Lifting / Handling for transport procedures.

TROUBLESHOOTING

CAUTION: Extreme caution must be exercised any time access is made into areas housing electrical components. Repair, replacement or adjustment of components in these areas must only be done by qualified technicians familiar with electrical circuitry and the operation of the chamber.

Always make a visual inspection of the incubator and control console when troubleshooting. Look for loose or disconnected wires or tubing, which may be the source of the trouble.

NOTE: IF BREAKERS TRIP REPEATEDLY, OR FUSES BLOW REPEATEDLY, CALL QUALIFIED SERVICE PERSONNEL.

TEMPERATURE

Temperature too high

- 1/ controller set too high-see section 5.3
- 2/ controller failed on – call Customer Service
- 3/ wiring error – call Customer Service

Display reads "HI" or "400"+

probe is unplugged, is broken or wire to sensor is broken – trace wire from display to probe; move wire and watch display to see intermittent problems

Chamber temp spikes over set point and then settles to set point

recalibrate – see section 5.4

Temperature too low

- 1/ OTP too low – see section 5.8.E
- 2/ controller set too low – see section 5.3
- 3/ unit not recovered from door opening – wait for display to stop changing
- 4/ unit not recovered from power failure or being turned off – incubators will need 24 hours to warm up and stabilize
- 5/ element failure – see if heating light is on; compare current draw to data plate
- 6/ controller failure – confirm with front panel lights that controller is calling for heat
- 7/ OTP failure – confirm with front panel lights that it is operating correctly
- 8/ wiring problem – check all functions and compare wiring to diagram in section 10.0, especially around any areas recently worked on
- 9/ loose connection – check shadow box for loose connections

Display reads "LO"

- 1/ sensor is plugged in backwards – reverse sensor wires to controller
- 2/ if ambient room temperature is lower than range of unit – compare set points and ambient temperature to rated specifications in section 9.0

Unit will not heat over a temperature that is below set point

- 1/ confirm that fan is moving and that amperage and voltage match data plate – check fan motor motion in shadow box and feel for air movement in chamber
- 2/ confirm that set point is set high enough ,turn OTP counterclockwise and see if heating light or safety light comes on
- 3/ check connections to sensor
- 4/ check calibration – using independent certified reference thermometer, follow instructions in sections 5.4 and 5.8

Unit will not heat up at all

- 1/ verify that controller is asking for heat by looking for controller light – if pilot light is not on continuously during initial start up there is a problem with the controller
- 2/ check amperage – amperage should be virtually at maximum rated (data plate) amperage
- 3/ do all controller functions work?
- 4/ is the OTP set high enough? – for diagnostics, should be fully clockwise with the pilot light never on
- 5/ has the fuse/circuit breaker blown?

Indicated chamber temperature unstable

- 1/ ± 0.1 may be normal
- 2/ is fan working? – remove top panel and verify movement of cooling fan in center of shadow box
- 3/ is ambient room temperature radically changing – either door opening or room airflow from heaters or air conditioning? – stabilize ambient conditions
- 4/ sensor miss-located, damaged, or wires may be damaged - check mounts for control and OTP sensors, then trace wires between sensors and controls
- 5/ calibration sensitivity – call Customer Service
- 6/ OTP set too low – be sure that OTP set point is more than 5 degrees over desired Main set point; check if OTP pilot is on continuously; turn controller knob completely clockwise to see if problem solved then follow instructions in section 5.8.E for correct setting.
- 7/ electrical noise – remove nearby sources of RFI including motors, arcing relays or radio transmitters
- 8/ bad connection on temperature sensor or faulty sensor – check connectors for continuity and mechanical soundness while watching display for erratic behavior; check sensor and wiring for mechanical damage
- 9/ bad connections or faulty capacitor – check connectors for mechanical soundness and look for corrosion around terminals or signs of arcing or other visible deterioration

Will not maintain set point

- 1/ assure that set point is at least 5 degrees over ambient room temperature
- 2/ see if ambient is fluctuating

Display and reference thermometer don't match

- 1/ calibration error – see section 5.4
- 2/ temperature sensor failure – evaluate if pilot light is operating correctly
- 3/ controller failure – evaluate if pilot light is operating correctly
- 4/ allow at least two hours to stabilize
- 5/ verify that reference thermometer is certified

Can't adjust set points or calibration

- 1/ turn entire unit off and on to reset
- 2/ if repeatedly happens, call Customer Service

Calibrated at one temperature, but not at another

This can be a normal condition when operating temperature varies widely. For maximum accuracy, calibration should be done at or as close to the set point temperature.

CO2 LEVEL

Overshoots set point but stabilizes - display and Fyrite match

- 1/ turn set point up and down to see if solenoid valve works by feeling and listening to valve
- 2/ recalibrate with Fyrite, see section 5.11 and section 7.0
- 3/ fan not operating correctly
 - a- fan motor stopped
 - b- fan blade fell off
 - c- wrong fan blade installed or mounted backwards
- 4/ wrong restrictor installed or missing altogether
- 5/ tank pressure too high, see section 5.9.A
- 6/ CO2 sensor partially plugged with dirt or condensation
- 7/ regulator set wrong, see section 5.9
- 8/ incubator too heavily loaded
- 9/ incubator being operated without shelving
- 10/ CO2 attached to sample port instead of "CO2 IN" fitting
- 12/ if display and Fyrite does not match, see Troubleshooting Co2-5

Overshoots set point and continues to rise - display and Fyrite match

- 1/ debris in solenoid causing it to leak continuously
- 2/ solenoid failed while open
- 3/ unit miss-wired
- 4/ controller output failed or shorted
- 5/ CO2 sensor or interface failure
- 6/ CO2 sensor plugged by debris or condensation

Rises very slowly

- 1/ restrictor partially plugged
- 2/ filter overly dirty or partially plugged
- 3/ CO2 tank regulator set too low, see section 5.9
- 4/ hose kinked or leaking
- 5/ poor door seal
- 6/ CO2 tank contains mixed gas, not 100% medical grade CO2

Never rises

- 1/ CO2 tank empty
- 2/ solenoid failed while closed
- 3/ CO2 controller output failed while open
- 4/ restrictor plugged
- 5/ CO2 hose blockage
- 6/ CO2 filter plugged
- 7/ set point is at 0.0 and has not been reset, see section 5.3 and 5.4

Display and Fyrite reading do not match

- 1/ calibration error - clear chamber for 12 hrs and confirm at "0"
- 2/ turn set point up and down to see if solenoid valve works by feeling and listening to valve
- 3/ assure that IR tube block is completely tight, 24 series
- 4/ if display is varying when doing Fyrite test, have a leak on IR base or mounting plate, or CO2 is injecting at time of test.
- 5/ take voltage reading across sensor leads 0=0V and 20=1V
- 6/ assure that sensor is properly mounted
- 7/ DISPLAY LOWER THAN FYRITE: air leak around the CO2 sensor or the fan motor shaft
- 8/ DISPLAY HIGHER THAN FYRITE: leak in tube from sample port to chamber, verify with liquid soap solution.
- 9/ DISPLAY HIGHER THAN FYRITE: Fyrite fluid worn out, replace fluid with Fyrite refill.
- 10/ DISPLAY HIGHER THAN FYRITE: poor Fyrite sampling technique, see section 6.0
- 11/ DISPLAY HIGHER THAN FYRITE: defective Fyrite pump or tubing
- 12/ CO2 sensor, interface or controller failure

Is unstable – display or actual reading varies around set point

- 1/ confirm that fan is working
- 2/ check for air leak around CO2 sensor or fan motor shaft
- 3/ magnetic field interference
- 4/ door not sealing properly
- 5/ atmospheric pressure fluctuations
- 6/ defective Fyrite pump or tubing
- 7/ CO2 entering chamber too quickly
 - a- tank pressure too high
 - b- regulator set wrong
 - c- restrictor incorrect or missing
- 8/ electronic problem with CO2 sensor, interface or controller
- 9/ top of unit exposed to cold air drafts
- 10/ unit being operated without shadow box cover
- 11/ electrical interference – atmospheric or conducted
- 12/ incubator too heavily loaded

Can't adjust set points or calibration - "locked up"

- 1/ turn unit off and on to reset processor in controller
- 2/ if repeatedly happens, call Customer Service

Feeding continuously or abnormally high CO2 usage

- 1/ do decay test: if more than 1% decay in an hour, check for leak : door gasket tightness, motor inlet to chamber, sensor or probe inlet to chamber, chamber access tube plugged
- 2/ check output signal from controller
- 3/ check solenoid valve for correct operation
- 4/ leak in plumbing including between regulator and CO2 tank
- 5/ door being opened too often

Won't hold calibration – Fyrite reading varies but display stable

- 1/ atmospheric pressure fluctuations
- 2/ top of unit exposed to cold air drafts
- 3/ unit being operated without shadow box cover in place
- 4/ condensation collecting on CO2 sensor
- 5/ CO2 sensor or interface failure
- 6/ unit incorrectly calibrated, see section 5.4
- 7/ taking Fyrite reading too soon after the door has been opened
- 8/ air leak around CO2 sensor mounting plate

MECHANICAL

Door not sealing

- 1/ check physical condition of gasket
- 2/ confirm that door latch pulls door in tightly
- 3/ assure that gasket is in original location

Motor doesn't move

- 1/ if shaft spins freely: check connections to motor and check voltage to motor;
- 2/ if shaft rubs or is frozen, relieve binding and retest

Motor makes noise

- 1) Make sure that the fan or blower wheel is not contacting its housing. Adjust the motor mounting bracket position to re-center the fan or blower wheel, if necessary.
- 2) Check the fan or blower wheel for damage or out of balance condition. Replace the fan or blower wheel if it is damaged or out of balance.
- 3) Turn the motor shaft to make sure that it spins freely. If it binds or the bearings make a rubbing or scrapping sound then replace the motor.

Solenoid valve buzzing

After removing solenoid clean with alcohol for carbon build-up then blow out. Check valve seat or channels for contamination. Check CO2 filter and/or grade of CO2 used. If seat is worn, replace.

OTHER

Controller on at all times - "locked-up"

- 1/ turn unit off and on to reset
- 2/ if cannot change any condition on the front panel, call Customer Service

Front panel displays are all off

- 1/ Check for wire damage.

Unit or wall fuse/circuit breaker is blown

- 1/ check wall power source
- 2/ compare current draw and compare to specs on data plate
- 3/ see what other loads are on the wall circuit

Unit will not turn on

- 1/ check wall power source
- 2/ check fuse/circuit breaker on unit or in wall
- 3/ see if unit is on, e.g., fan or heater, and just controller is off
- 4/ check all wiring connections, esp. around the on/off switch

Unit is smoking – Out of box

Put unit under vent and run at full power for one hour.

Contamination in chamber

- 1/ see cleaning procedure in Installation section 3.0
- 2/ develop and follow Standard operating procedure for specific application; include definition of cleaning technique and maintenance schedule.

NOTE: THE TEMPERATURE OF THE GLASS DOOR AND THE INTERIOR CHAMBER WALLS MUST BE EQUAL TO PREVENT CONDENSATION.

The incubator is designed so that no internal electrical servicing should be required under normal conditions. If electrical servicing is necessary, it should be performed by qualified service personnel. **FOR PERSONAL SAFETY, ALWAYS DISCONNECT THE POWER BEFORE SERVICING.**

PARTS LIST

Description	115V	220V
Blower Wheel	2600519	2600519
CO2 Display Board	1750559	1750735
CO2 Filter	100199	100199
CO2 Pre-Heater	100909	103328
CO2 Sensor w/ Interface	9700501	9700501
Cord Set	1800510	104192
Cord Set - European	NA	1800500
Data Logger	1750537	1750737
Door Heater	103068	103068
Door Switch	X1000022	X1000022
Element Coil	2350503	120071
Fan Blade	103130	103130
Fan Motor	4880508	4880507
Glass Door Latch	9530516	9530516
Gasket, 8 feet	3450648	3450648
HEPA Filter	2800517	2800517
Humidity Display	1750736	1750736
Humidity Pan	5120587	5120587
Humidity Sensor	9900522	9900523
IR Transmitter	8320510	8320510
Leveling Feet	200129	200129
Limit Thermostat	103557	103557
Overtemperature Safety	100001	100001
Pod Control 4 Loop	1750725	1750725
Power Inlet Fuse	103558	103555
Power Inlet	104179	104179
Power Switch	7850532	7850532
Shelf	5121109	5121109
Solenoid Valve	8600528	8600529
Temperature Display Board	1750557	1750734
Water Level Switch	210004	210004

UNIT SPECIFICATIONS

Weight	Shipping	Net
3507 & 3517	280 lbs.	221 lbs.
3514 & 3524	560 lbs.	442 lbs.


Dimensions	Exterior WxDxH (IN.)	Interior WxDxH (IN.)
3507 & 3517	26.5 x 26 x 40	20.5x20x28
3514 & 3524	26.5 x 26 x 80	20.5 x 20x28 each chamber

Capacity	Cubic Feet
3507 & 3517	6.7
3514 & 3524	6.7 each chamber

Temperature	Range	Uniformity	CO2
3507 & 3517	Amb. 8°C to 60°C	±0.25°C @ 37°C	0 – 20%
3514 & 3524	Amb. 8°C to 60°C	±0.25°C @ 37°C	0 – 20%


Model	Amp	Voltage	Hertz
3507	5 amp	120V	60 Hz
3507-2	3 amp	220V	50/60 Hz
3517	5 amp	120V	60 Hz
3517-2	3 amp	220V	50/60 Hz
3514	5 amp	120V	60 Hz
3514-2	3 amp	220V	50/60 Hz
3524	5 amp	120V	60 Hz
3524-2	3 amp	220V	50/60 Hz

FIGURE 5
WIRING DIAGRAM
3507 & 3514
(TC Sensor)

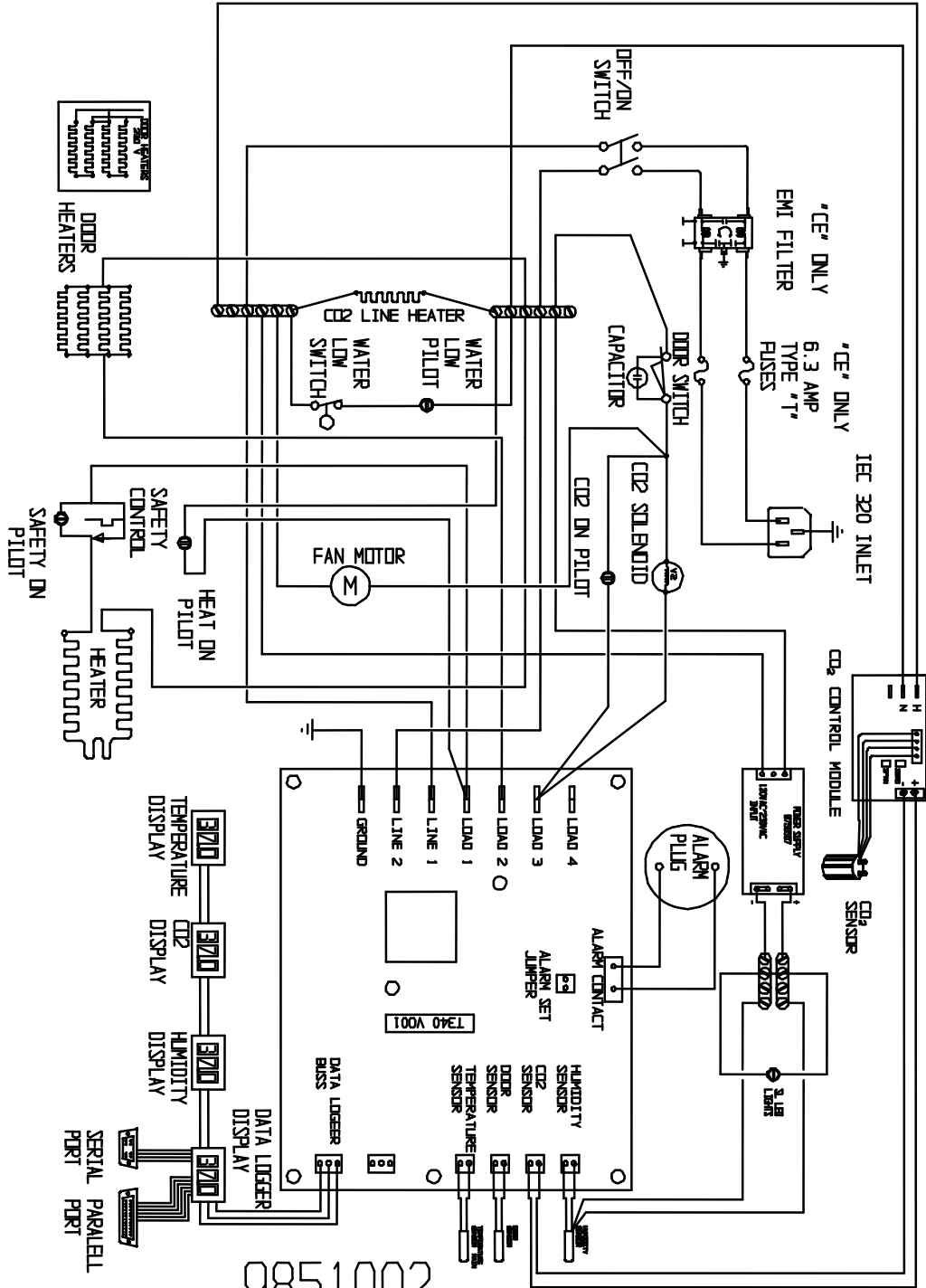

9851002

**FIGURE 6
WIRING DIAGRAM
3517 & 3524
(IR Sensor)**

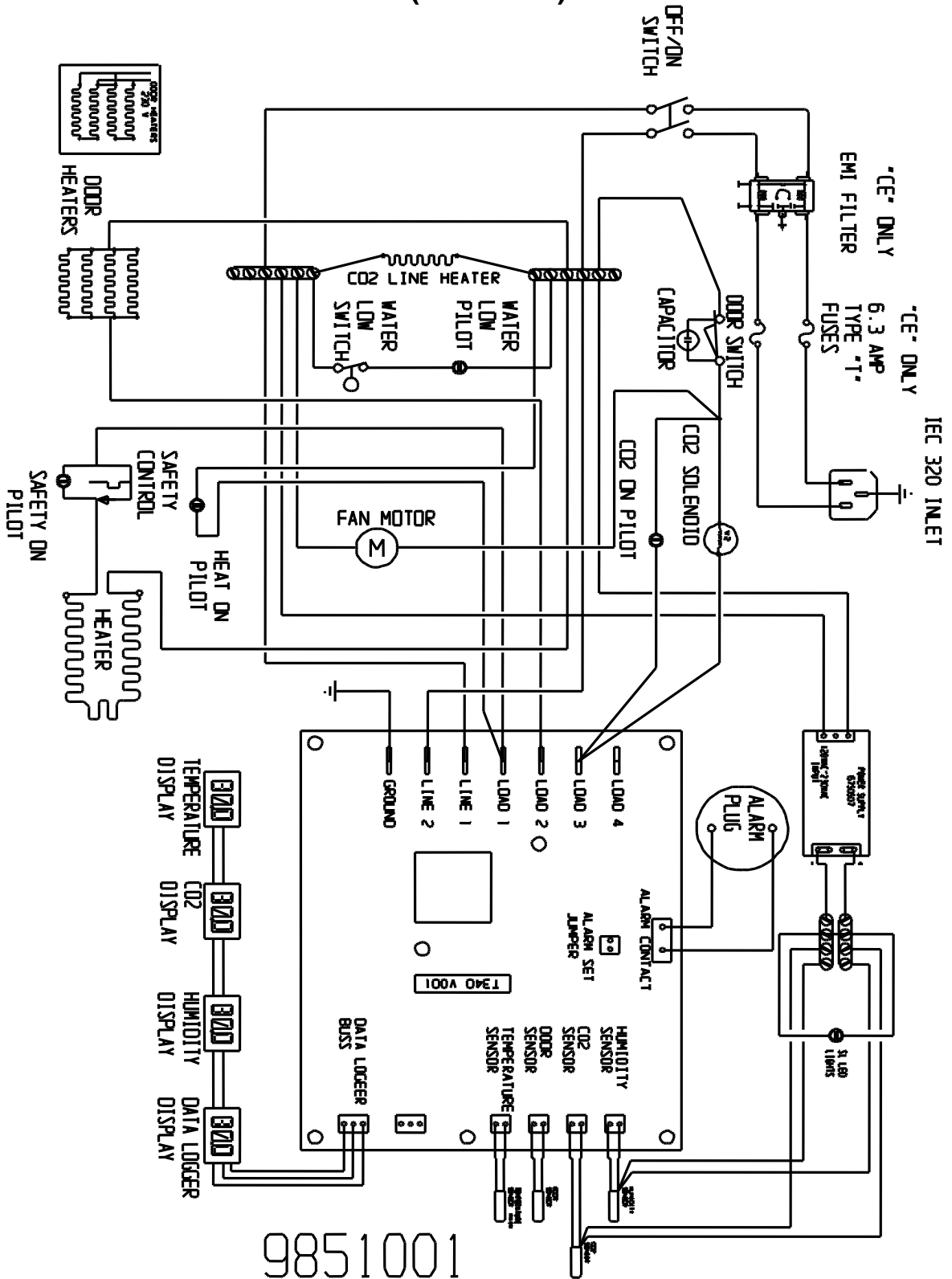

9851001